

Spinsolve® 60

60 MHz benchtop NMR

The Magritek Spinsolve 60 NMR spectrometer sets a new standard for 60MHz benchtop NMR instruments. The Spinsolve 60 spectrometer provides high quality ^1H and ^{19}F NMR spectra in seconds, right where it is needed - at your laboratory bench.





Features

- ^1H and ^{19}F nuclei
- 1D and 2D NMR experiments
- Compact, benchtop size and weight
- No spinning or compressed gas
- No cryogenes

Applications

- Research laboratories
- Industrial QA/QC laboratories
- Undergraduate education

Example spectra from the Spinsolve 60 benchtop NMR

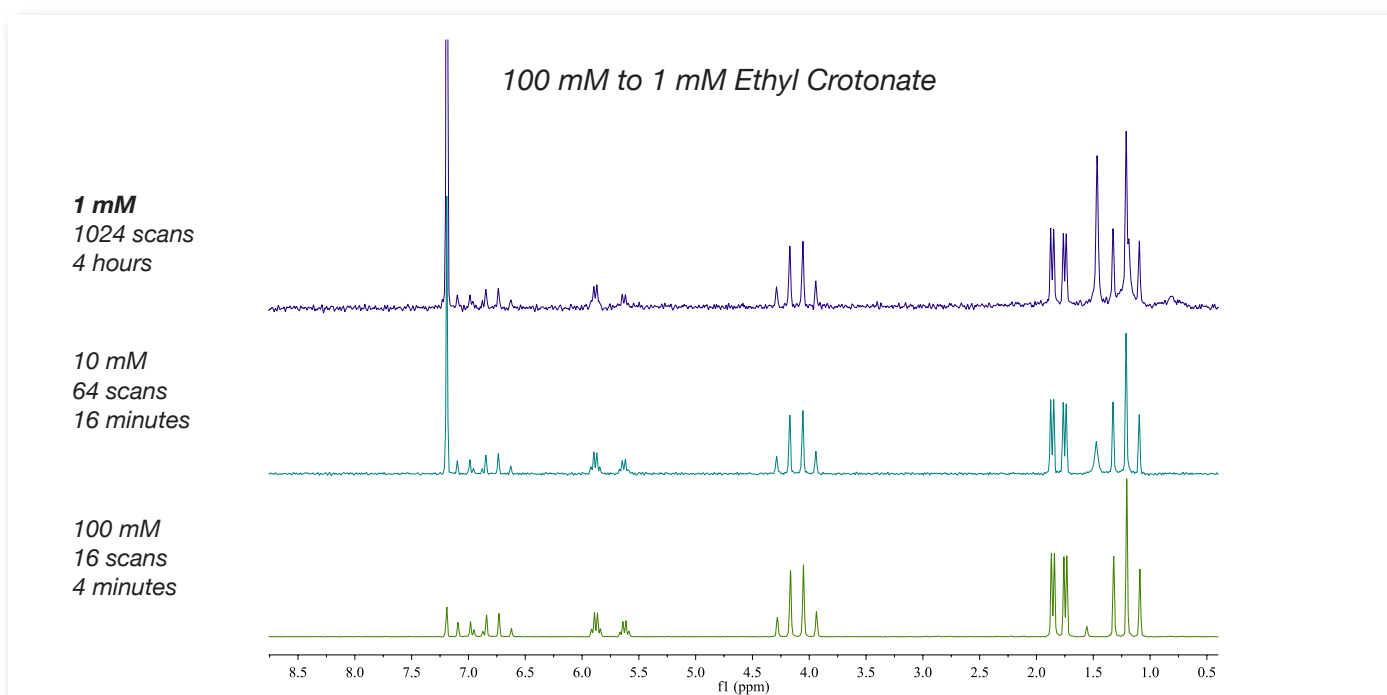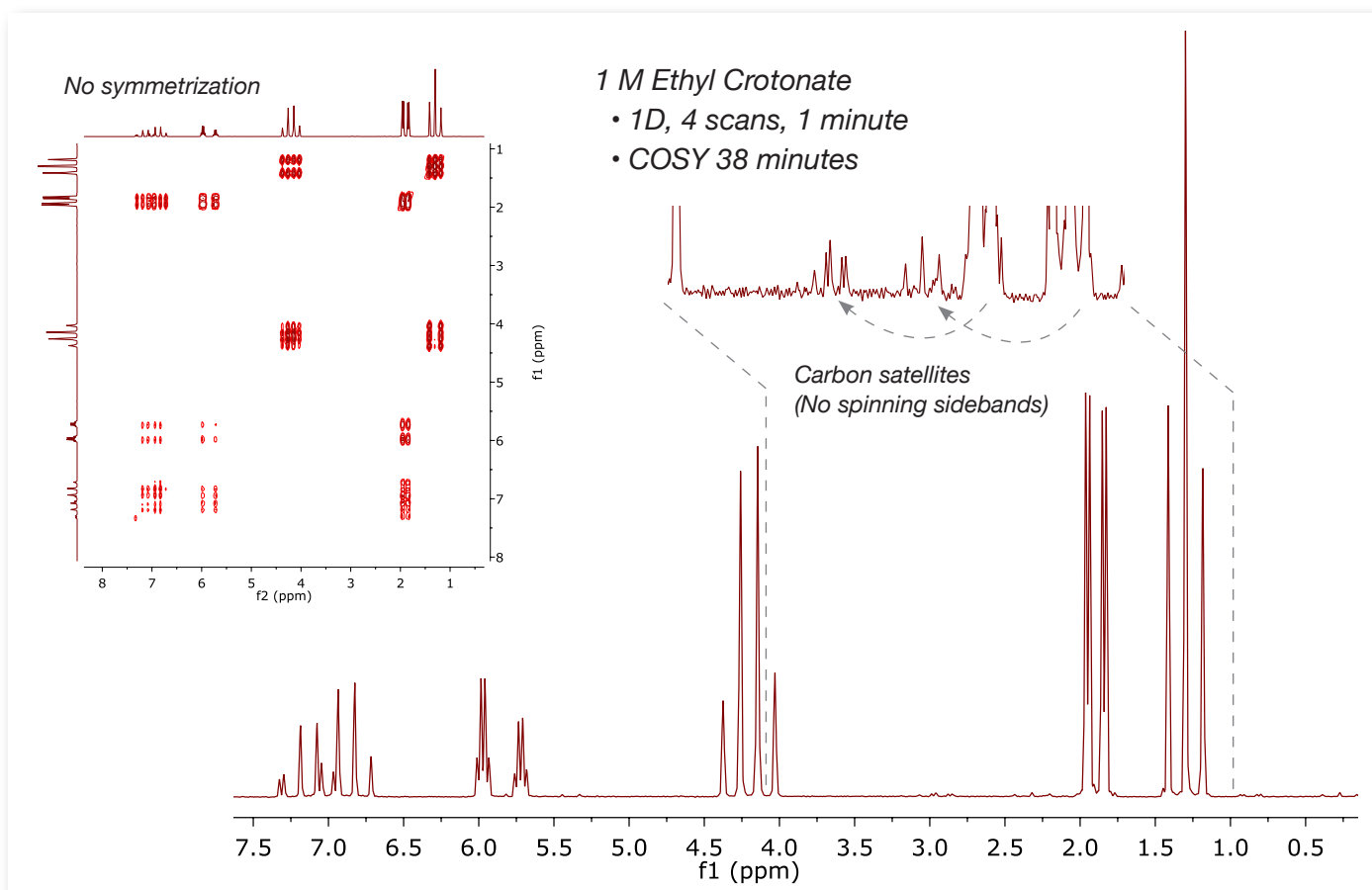

Software


The Spinsolve software is beautifully simple and easy to use, with a clean and intuitive user interface.

Easy experiment selection in the upper menu

Single button start

Progress bar

One-click processing buttons

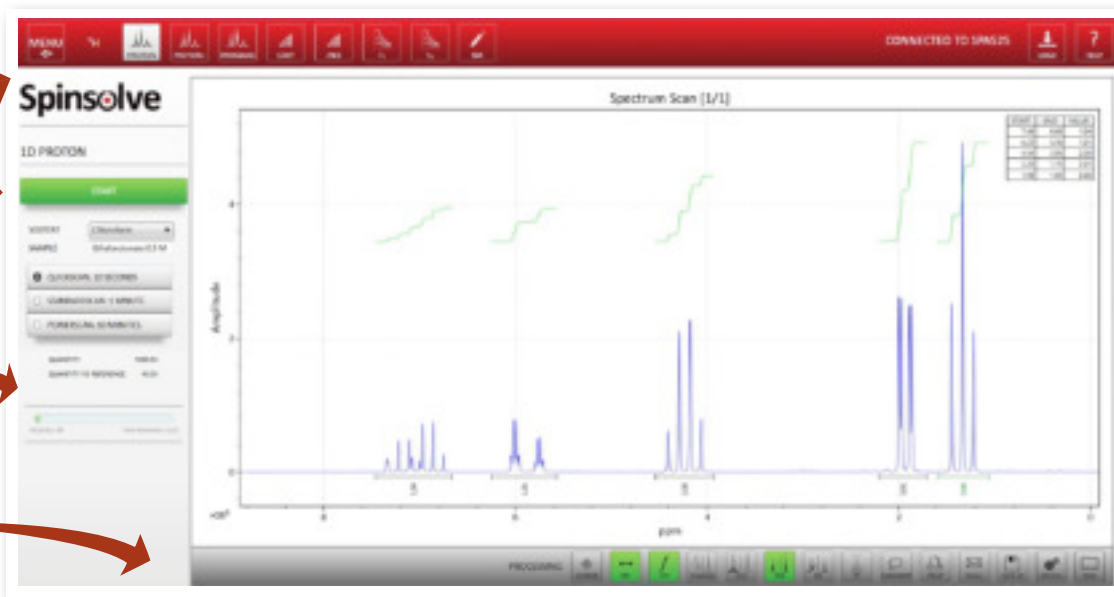

Simple menu structure


1. Click to choose nucleus
2. Click to choose experiment
3. Click Start (watch status on the progress bar)
4. Click any processing you wish to apply



Automate experiments with scripts

Using scripts the user can queue a series of experiments to run sequentially. A range of script templates are provided for easy modification. Scripts are displayed graphically to provide a clear picture of the sequence.

This script shown here performs a 1D Proton experiment, followed by a JRES and then a COSY.

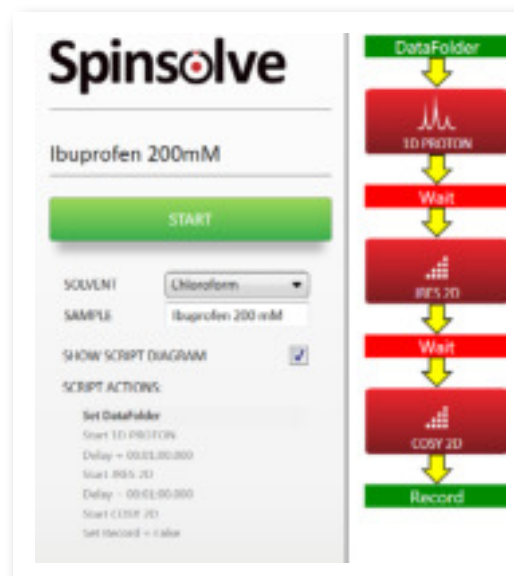


Spinsolve 60



Specifications

- Nuclei: ^1H , ^{19}F
- Operating frequency: 60 MHz (^1H)
- ^1H 50% Linewidth: < 0.5 Hz
- ^1H 0.55% Linewidth: < 20 Hz
- ^1H Sensitivity: $>180:1$ for 1% Ethyl Benzene
- Operating Temperature Range: 20°C to 25°C (68°F to 77°F)
- Dimensions: 58 x 43 x 40 cm (23" x 17" x 16")
- Weight: 60 kg (120 lb)
- Stray Field: < 2 G all around system
- Power requirement: 110-240V AC

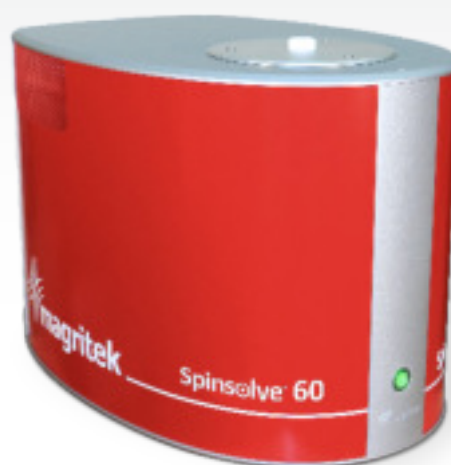

Pulse sequences available on the Spinsolve 60 spectrometer

Proton	Fluorine
1D	1D
Paramagnetic	Paramagnetic
2D COSY	2D F - COSY
2D TOCSY	2D F - JRES
2D JRES	2D FH - COSY
T_1 , T_2	Reaction Monitoring
Reaction Monitoring	

Contact us now for a quote, to request a demo or to measure your samples

Email: sales@magritek.com

Website: www.magritek.com/contact-us

GERMANY +49 241 9278 7270

UNITED KINGDOM +44 7468 529 615

UNITED STATES +1 855 667 6835

NEW ZEALAND +64 4 477 7096

For a complete listing of our global offices and distributors visit: www.magritek.com/about-us/distributors/