## Artemisinin

# Spinsolve 90

Artemisinin is a widely used drug in the standard treatment of malaria. It is extracted from the plant *Artemisia annua*, sweet wormwood, but can also be produced in a semi-synthetic fashion. Figure 1 shows the <sup>1</sup>H NMR spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> measured in a single scan taking 10 seconds to acquire.

#### 1D Proton spectrum

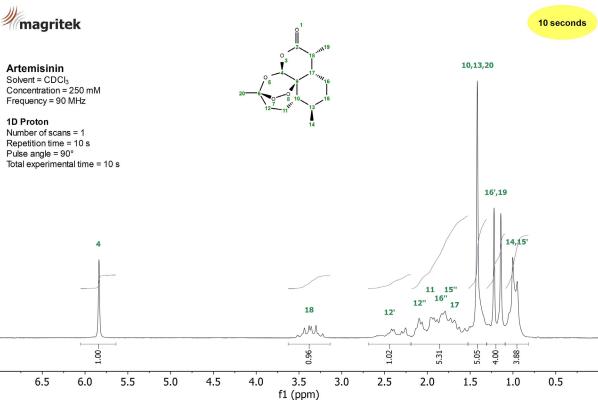


Figure 1: <sup>1</sup>H NMR spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> measured on a Spinsolve 90 MHz system in a single scan.

### 1D Carbon spectrum

Figure 2 shows the <sup>13</sup>C NMR spectrum of 250 mM Artemisinin in CDCl<sub>3</sub> acquired using NOE polarization transfer from <sup>1</sup>H to <sup>13</sup>C and <sup>1</sup>H decoupling. The 1D Carbon experiment using NOE is sensitive to all <sup>13</sup>C nuclei in the sample. It clearly resolves all the expected resonances.

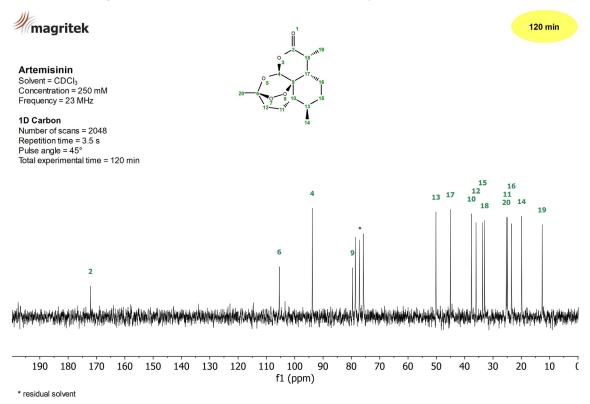


Figure 2: <sup>13</sup>C NMR spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> measured on a Spinsolve 90 MHz system in 120 minutes.



### 2D COSY spectrum

# Spinsolve 90

magritek

The 2D COSY experiment allows one to identify coupled <sup>1</sup>H nuclei as they generate cross peaks out of the diagonal of the 2D data set. In Figure 3 a large number of cross peaks can be nicely observed. For example, the protons at position 4 and 17 (dark blue) couple with each other. Furthermore, proton 18 couples with proton 17 (cyan) and 19 (pink).

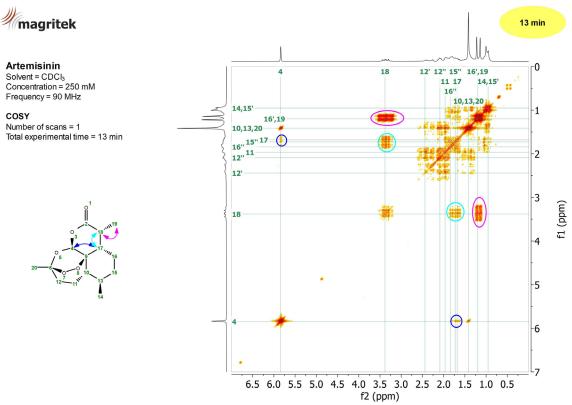


Figure 3: <sup>1</sup>H 2D COSY experiment of a 250 mM Artemisinin sample in CDCl<sub>3</sub> acquired in 13 minutes on a Spinsolve 90 MHz system.

#### 2D HSQC-ME

The HSQC is a powerful sequence widely used to correlate <sup>1</sup>H with the one-bond coupled <sup>13</sup>C nuclei. The Spinsolve is equipped with a multiplicity edited version (HSQC-ME) of this method. It provides the editing power of the DEPT-135 sequence, which is useful to differentiate the signals of CH<sub>2</sub> groups (blue) from CH and CH<sub>3</sub> groups (red). Figure 4 shows the HSQC-ME spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> acquired in 8 minutes. The measurement time was optimized applying NUS (non uniform sampling).

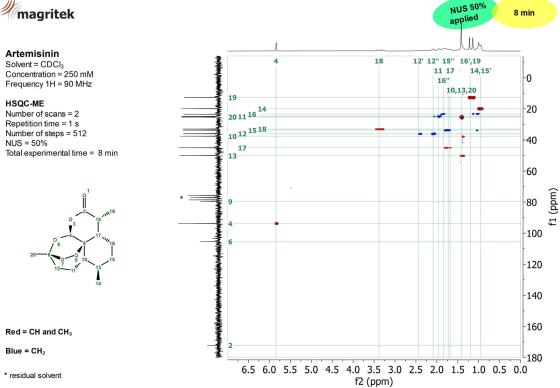


Figure 4: HSQC-ME spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> showing the correlation between the <sup>1</sup>H (horizontal) and <sup>13</sup>C (vertical) signals.

### 2D HMBC

# Spinsolve 90

To obtain long-range <sup>1</sup>H-<sup>13</sup>C correlations through two or three bond couplings, the Heteronuclear Multiple Bond Correlation (HMBC) experiment can be used. Figure 5 shows the HMBC spectrum of a 250 mM Artemisinin sample measured in 34 minutes on our Spinsolve 90 MHz. As an example, the long-range correlation of protons 19 with carbons 2, 17 and 18 are marked. The experiment shows the correlation with quaternary carbons, too.

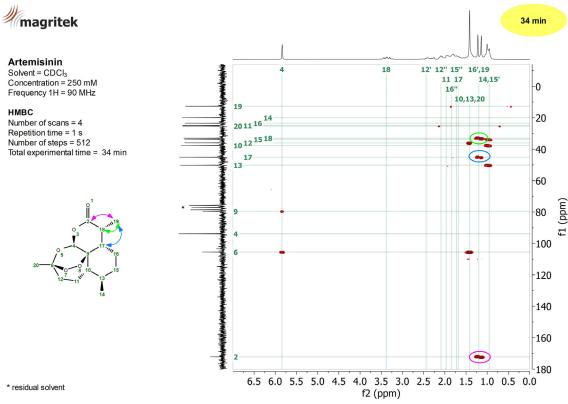


Figure 5: HMBC spectrum of a 250 mM Artemisinin sample in CDCl<sub>3</sub> showing the long-range couplings between <sup>1</sup>H and <sup>13</sup>C nuclei.

